

Cat.# R1003

1-Shot Digital ECL HRP Substrate Solution

Kit Component:

1-Shot Digital ECL HRP Substrate Solution (125 mL)

Instruction

Storage: 4°C

1. Perform Western blotting according to users' normal protocols.
2. Rinse the membrane with dI water for 30 seconds.
3. Apply the solution directly on the membrane and cover the entire surface;
4. Incubate at room temperature for 2 minutes, and then drain the surface remaining solution, and lay face-up on *KwikQuant* Imager's sample tray.
5. Acquire the blot images using *KwikQuant* Imager.