

Clean Western Buffer (Cat.# R2001)

Blocking membrane is not needed if this buffer used to detect bands on the membrane. Please see product # R1100's protocol.

Clean Western Buffer produces cleaner blots than regular TBST does.

Use this buffer to replace TBST in blocking, primary antibody blotting, secondary antibody blotting, and every washing step.

IMPORTANT: resins the blot with dI-water before applying ECL substrate solution. This buffer contains a compound that inhibits HRP.

Preparation: Add component A and B in 3.5L de-ionized water. Shake 10-15 minutes to dissolve all components. Store at room temperature.